

OLYMPUS®

Olympus is about life. About photographic innovations that capture precious moments of life. About advanced medical technology that saves lives. About information- and industry-related products that make possible a better living. About adding to the richness and quality of life for everyone. Olympus. Quality products with a **FOCUS ON LIFE**

POLARIZING MICROSCOPE

BX51-P


BX2 SERIES

UIS
UNIVERSAL
INFINITY SYSTEM




Unsurpassed optics render polarized light images sharper than ever before.

Olympus is proud to introduce the BX51-P, the new polarizing microscope with superb performance in polarized light. It's a breakthrough combination of world-renowned UIS infinity-corrected optics and Olympus original optical-design. Olympus has also devised a new, extended line of compensators to make the BX51-P versatile enough to handle observation and measuring applications in virtually any field.

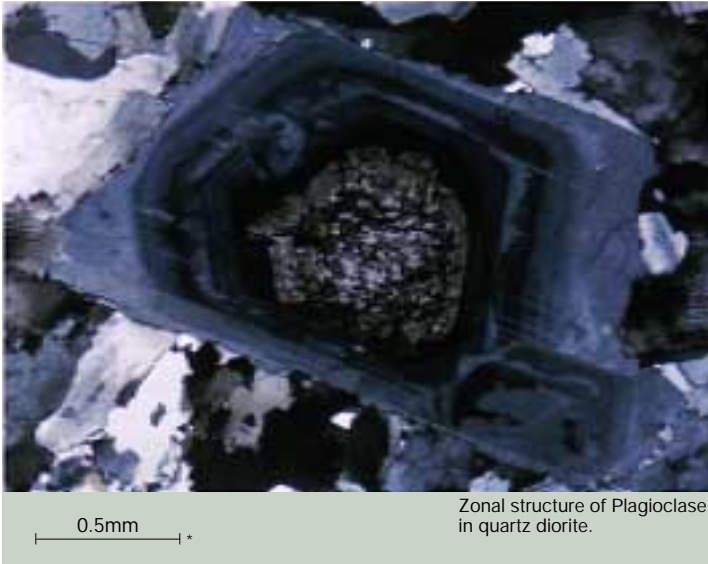

Zonal structure of Plagioclase in quartz diorite.

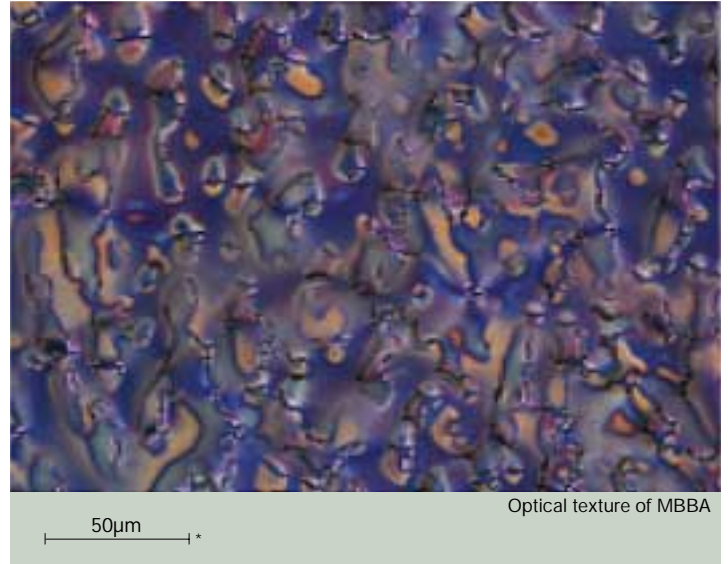

Optical texture of MBBA

**Scales indicate actual size of samples*

UIS provides outstanding system expandability.

By maximizing the advantages of infinity correction, the UIS optical system prevents the deterioration of optical microscope performance and eliminates magnification factors, even when polarizing elements like analyzers, tint plates or compensators are introduced into the light path. Maintaining the highest level of system flexibility, the BX51-P also accepts intermediate attachments available for the BX2 Series system microscopes, as well as video cameras and photomicrographic systems.



BX51-33P
(Conoscopic/Orthoscopic version)



BX51-33PO
(Orthoscopic version)




Enhanced operational ease and outstanding optical performance.

- With the U-CPA conoscopic observation attachment, the changeover between orthoscopic and conoscopic observation methods is simple and quick — just slide the Bertrand lens control knob in or out.
- The circular rotatable stage has two centering knobs and allows smooth sample rotation. A settable click stop every 45 degrees enables accurate observation and measurement.
- An ergonomic Y-shaped frame ensures comfortable, efficient observation with less fatigue, even over prolonged periods.
- Standard field number F.N. 22 represents a 21% increase in field size, covered with the conventional F.N. 20.
- The high intensity 12V, 100W halogen light source combined with the UIS optical system and its high transmittance result in exceptionally bright and sharp images.
- Mounting an attachable cross-movement mechanical stage (U-FMP) onto the circular rotatable stage makes for improved observation efficiency. Interference between the mechanical stage and the objectives is eliminated, so that images of superb quality can be effortlessly observed with all objective magnification.




U-FMP

BX51-P system diagram

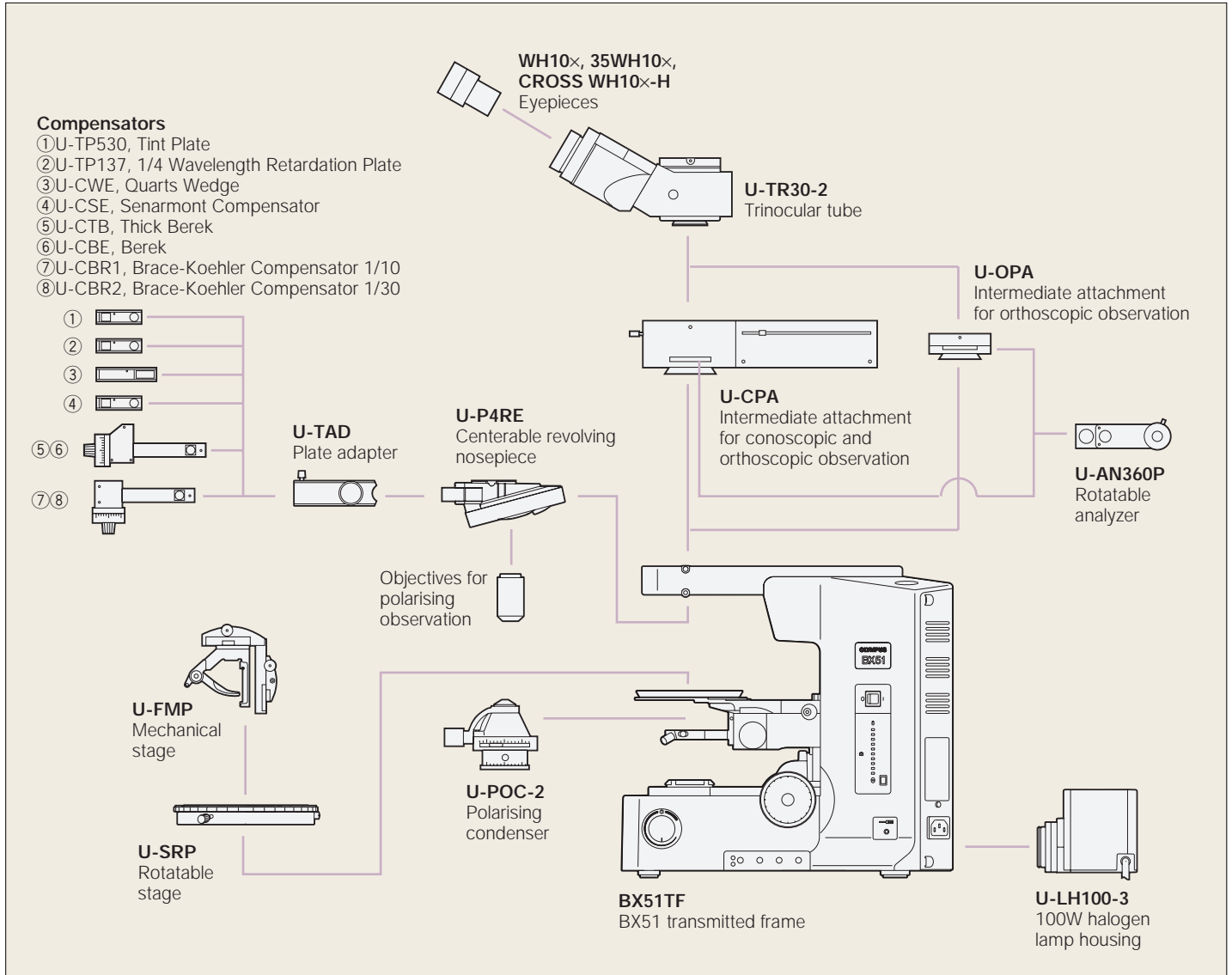

BX51-P dimensions

(unit: mm)

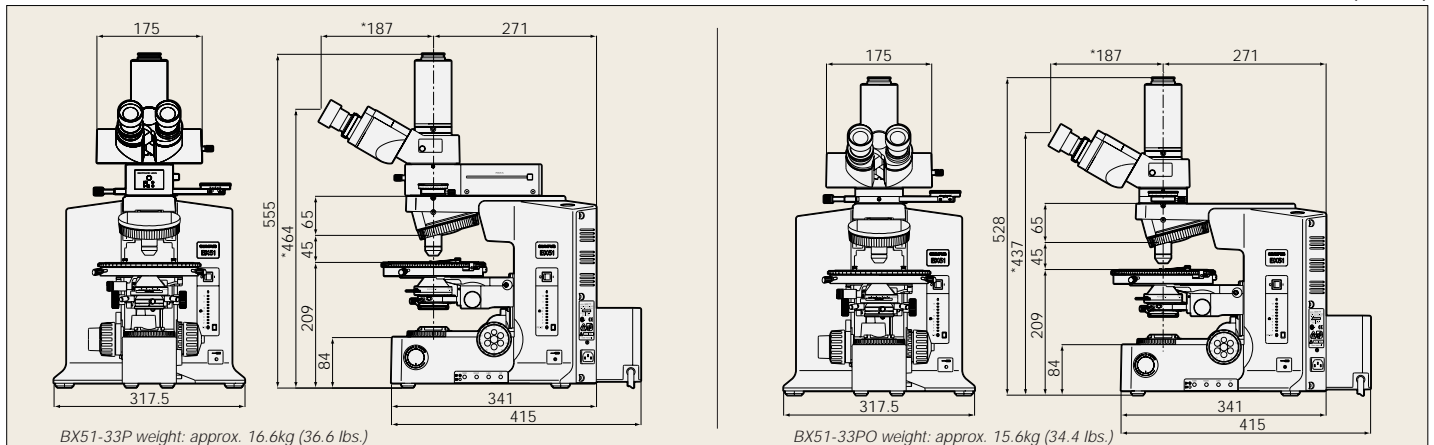

Power consumption: 250VA Note: The length marked with an asterisk (*) may vary according to the interpupillary distance setting. Distance for figure is 62mm.

An upgrade in polarization characteristics.

Thanks to Olympus' sophisticated design and manufacturing technology, new ACH-P and UPLFL-P strain-free objectives reduce internal strain to an absolute minimum. Olympus has also totally redesigned its polarizers and polarizing condensers to further enhance performance in polarized light. This means a higher EF* value, resulting in unmatched image contrast. To meet a diversity of research and application requirements, the universal UPLFL-P Series objectives have been designed to accommodate a wide range of observation methods, including Nomarski DIC and fluorescence microscopy, in addition to polarized light observation.

*The EF (extinction factor) is the brightness ratio between parallel and crossed pol-filters. The higher the EF value, the better the extinction.



UPLFL-P Series



ACH-P Series

UPLFL-P Series

Item		N.A.	W.D.
UPlan FL	4xP	0.13	13.0mm
UPlan FL	10xP	0.30	3.1mm
UPlan FL	20xP	0.50	1.6mm
UPlan FL	40xP	0.75	0.51mm
UPlan FL	100xOP	1.30	0.10mm

PL-P Series

Item		N.A.	W.D.
Plan FL	4xP	0.1	22.0mm

ACH-P Series

Item		N.A.	W.D.
ACH	10xP	0.25	6.1mm
ACH	20xP	0.40	3.0mm
ACH	40xP	0.65	0.45mm
ACH	100xOP	1.25	0.13mm



①U-TP530 ②U-TP137 ③U-TAD ④U-CBR1 ⑤U-CBR2 ⑥U-CWE ⑦U-CSE ⑧U-CBE ⑨U-CTB

An extensive range of compensators is available.

Six different compensators are available for the BX51-P microscope, allowing measurement of various retardation levels, ranging from 0 to 20λ. For easier

measurement, the direct readout method is featured. Higher image contrast can be attained by using a Senarmont* or Brace-Koehler compensator to change the retardation level in the entire field of view.

* Used with monochromatic green filter, IF546 or IF550.

Measuring range of compensators

Compensator	Measurement range	Applications
Thick Berek (U-CTB)	0-11,000nm (20λ)	Measurement of high retardation level (R' > 3λ), (crystals, macromolecules, fiber, etc.)
Berek (U-CBE)	0-1,640nm (3λ)	Measurement of retardation level (crystals, macromolecules, living organisms, etc.)
Senarmont compensator (U-CSE)	0-546nm (1λ)	Measurement of retardation level (crystals, living organisms, etc.) Enhancement of image contrast (living organisms, etc.)
Brace-Koehler compensator 1/10λ (U-CBR1)	0-55nm (1/10λ)	Measurement of low retardation level (living organisms, etc.)
Brace-Koehler compensator 1/30λ (U-CBE2)	0-20nm (1/30λ)	Enhancement of image contrast (living organisms, etc.)
Quartz wedge (U-CWE)	500-2,200nm (4λ)	Approximate measurement of retardation level (crystal, macromolecules, etc.)

*R= retardation level

For more accurate measurement, it is recommended that compensators (except U-CWE) be used together with the interference filter 45IF546.



①U-CPA ②U-P4RE ③U-AN360P ④U-OPA ⑤U-POC-2

Unmatched sharpness in orthoscopic and conoscopic observations.

With a U-CPA conoscopic observation attachment, changeover between orthoscopic and conoscopic observation is simple and quick.

Focusing of conoscopic images is easy and accurate. Employing a Bertrand field stop makes it possible to obtain consistently sharp and clear conoscopic images.

BX51-P specifications

Item	Conoscopic and orthoscopic observation (U-CPA)	Orthoscopic observation (U-OPA)
Polarized light intermediate attachment (U-CPA or U-OPA)	F.N.	22
	Bertrand lens	Focusable
	Bertrand field stop	ø3.4mm diameter (fixed)
	Engage or disengage Bertrand lens changeover between orthoscopic and conoscopic observation	Position of slider ● in Position of slider ○ out
	Analyzer slot	Rotatable analyzer with slot (U-AN360P)
Analyzer (U-AN360P)	360° dial-rotatable Rotatable minimum angle 0.1°	
Revolving centerable nosepiece (U-P4RE)	Quadruple, centerable attachable components: 1/4 wavelength retardation plate (U-TAD), tint plate (U-TP530) and various compensators can be attached using plate adapter (U-TAD)	
Stage (U-SRP)	Polarizing rotatable stage with 3-point centering function. 360° rotatable, lockable in any position, 360° graduated in 1° increments (minimum retardation resolution 6', using vernier scale) 45° click stop function ● Slide holder (U-SCB2) and mechanical stage (U-FMP) can be attached	
Condenser (U-POC-2)	Achromat strain-free condenser (U-POC-2), 360° rotatable polarizer with swing-out achromatic top-lens, Click stop at position "0°" is adjustable. N.A. 0.9 (top-lens in) N.A. 0.18 (top-lens out) Aperture iris diaphragm: adjustable from 2mm to 21mm diameters	



Specifications are subject to change without any obligation on the part of the manufacturer.

Olympus business areas

	Medical and health-care area
	Imaging and information area
	Industrial applications area



OLYMPUS OPTICAL CO., LTD.
 San-Ei building, 22-2, Nishi Shinjuku 1-chome, Shinjuku-ku, Tokyo, Japan
OLYMPUS OPTICAL CO. (EUROPA) GMBH.
 Postfach 10 49 08, 20034, Hamburg, Germany
OLYMPUS AMERICA INC.
 2 Corporate Center Drive, Melville, NY 11747-3157, U.S.A.
OLYMPUS SINGAPORE PTE LTD.
 491B River Valley Road, #12-01/04 Valley Point Office Tower, Singapore 248373
OLYMPUS OPTICAL CO. (U.K.) LTD.
 2-8 Honduras Street, London EC1Y 0TX, United Kingdom.
OLYMPUS AUSTRALIA PTY. LTD.
 104 Ferntree Gully Road, Oakleigh, Victoria, 3166, Australia